



DISS 6 Hardware

Digital image scanning system for scanning electron microscopes (SEM)

User manual

Document version: 2.1

Date of issue: 2026-03-04

Document language: EN

Translation of German original

Use Anybody working with/on the device has to have read and understood the relevant parts of this document.

Access The staff working with the device has to have constant access to this document to prevent handling errors and guarantee trouble-free operation.

Reproduction All information in this document is protected by copyright. This document must not be copied, distributed or changed in any way by anyone including business units or departments of point electronic GmbH without prior written approval by point electronic GmbH. Reproduction and use of this document are confined to internal purposes of the operator only.

Up-to-datedness Every endeavor has been made to make sure that the information contained in this document is complete and correct at the printing date. This document describes all units and functions known of at the current point of time.

Brand, company, or product names The brand, company, or product names used in this document are the brand, company, or product names of the respective manufacturers and/or owners.

Contact us point electronic GmbH
Erich-Neuß-Weg 15
06120 Halle (Saale)
Germany
Phone: +49 345 1201190
Email: info@pointelectronic.de
Web: <https://www.pointelectronic.de>

Product information Product: Digital image scanning system, 6th generation
Description: Digital image scanning system for scanning electron microscopes (SEM)
Name: PE-DISS6
Manufacturer: point electronic GmbH

Table of contents

1	Introduction	5
	User groups of this document	6
	Structure of this document	7
	Representations in this document	8
	Identification of the warning notices	9
	Topics relevant to service	10
2	Safety regulations	11
	Introduction	12
	Basic hazards	13
	Staff and qualifications	14
	Responsibility of the operating company	15
	Working in a safety-conscious manner	16
	Modifications and alterations	17
	Maintenance work	18
	Cleaning	20
	Environmental protection	21
3	System overview	22
	Intended use	23
	Improper use	24
	Scope of delivery	25
	Optional equipment	26
	System environment requirements	27
	Front panel	28
	Back panel	29
	Touch display	31
	Device labeling	34
4	Pinouts of the inputs and outputs	36
	Analog I/O	37
	Digital in	39
	Digital I/O	40
	AUX B1-4	41
	Mics M1-4	42
5	Installation and configuration	43
	Configuring the inputs and outputs	44
	Positioning	47
	Connecting to the power supply	48
	Installing	49
	Configuring TCP	51
6	Maintenance	54
	Replacing the fuses	55
7	Disposal	57
	Recycling and taking back of used equipment	58

8	Specifications	59
	Operating conditions	60
	Device specifications	61



1 Introduction

Chapter overview

Purpose This chapter contains information to simplify working with this document.

Contents This chapter contains the following information:

- › User groups of this document 6
- › Structure of this document 7
- › Representations in this document 8
- › Identification of the warning notices 9
- › Topics relevant to service 10

User groups of this document

Operators "Operators" are persons with right of disposal over the device using this right for purchase or other purposes.

Technical staff "Technical staff" are those persons that the operators entrust with tasks related to use and operation. Technical staff are trained by the operators to carry out their assigned tasks and are informed on the potential hazards that may arise due to improper handling.

Technical staff need to undergo a training by point electronic GmbH or a partner company of point electronic GmbH covering the following topics:

- Handling of the connected SEM
- Areas of application and operation of the device and the associated software
- Handling the connected peripherals
- Basic maintenance tasks and troubleshooting

Service technicians "Service technicians" are staff members whose specialist training, knowledge and experience as well as familiarity with the relevant safety regulations mean they are able to assess the tasks they have been entrusted with and anticipate potential hazards.

Service technicians may be:

- Employees of point electronic GmbH or
- Employees of partner companies of point electronic GmbH

Structure of this document

- Composition** This document is composed of chapters which are organized by technical aspects.
- Numbering** The chapters are numbered with Arabic numerals. Chapters may be composed of sections. Sections are numbered as second numbering level (e. g. 3.1). Sections are used to structure large chapters into sub-chapters.
- All pages of this document are numbered consecutively.
- Overviews** Each chapter and section contains an overview detailing contents and page numbers. This allows for direct access to a topic and independent use of single parts of this document.
- Related information** Related information within this document is marked by the note "Continuation next page ..." resp. "... continuation:". Please pay attention to the completeness of the information when copying parts of this document.
- Cross references** The content of this document is structured by topics. If further information on one topic may be found elsewhere in the user guide, the relevant chapter and page are pointed out.

Representations in this document

Illustrations Illustrations used in this document do not always contain all details or special cases. They only represent the relevant information.

Icons The following icons are used to mark certain information visually:

Icon	Meaning
	Notice For example:  The selected parameter will not be inserted in the parameter order.
	Reference to another part of this guidebook For example:  see "Cross references" on page 7
	Use of a tool For example:  Screwdriver TX 10

Notes Important notes are marked as follows:

NOTICE

Mind the notices in this document!

Notes explain relations that even for expert users might not be evident at first glance.

The neglect of a note is no direct security risk. However, it can lead to disturbances in the operating procedure.

Identification of the warning notices

Purpose This document includes warning notices that may lead to serious consequences if disregarded. Warning notices are not only listed in the “Safety regulations” chapter, but especially in places where hazards for people, equipment and operation may arise.

Identification of warnings There are three classes of warnings. These classes are indicated by specific signal words and colors. They include the following:

Signal word	Meaning
 DANGER	Warning notice, which if disregarded will probably or very likely result in death or serious injury.
 WARNING	Warning notice, which if disregarded may result in serious injury, permanent damage to health or serious property loss.
 CAUTION	Warning notice, which if disregarded may result in injury or property loss including financial losses due to operational impairment.
 ATTENTION	Warning notice, which if disregarded may result in property loss including financial losses due to operational impairment.

Topics relevant to service

Service contact For detailed information and questions about topics such as installation and configuration, please contact our service department at:

service@pointelectronic.de

Further documents Additional documents with descriptions and instructions for installation, configuration, and additional software tools are available in protected public areas.

If you require access to further documents, please contact our service department at:

service@pointelectronic.de

2 Safety regulations

Chapter overview

Purpose This chapter contains safety information relating to the protection of persons as well as safe and fault-free operation. All user groups of the hardware and software must be aware of and follow these safety provisions.

Contents This chapter contains the following information:

› Introduction	12
› Basic hazards	13
› Staff and qualifications	14
› Responsibility of the operating company	15
› Working in a safety-conscious manner	16
› Modifications and alterations	17
› Maintenance work	18
› Cleaning	20
› Environmental protection	21

Introduction

Reliable and safe operation The reliable and safe operation of the device depends on carefully implementing the operating, setup and maintenance tasks.

Observance of the safety instructions Always observe and follow the safety instructions and accompanying code of conduct when handling the device and the software. Always point out the instructions and code of conduct to staff working with the hardware and software.

Generally applicable safety regulations (such as accident prevention and environmental protection regulations, etc.) must also be observed.

Consequences arising from failure to observe the safety instructions Failure to observe the safety instructions can result in death or serious injury of personnel and damage or destruction of the hardware components.

Basic hazards

Definition The device meets the current state of the art as well as the applicable safety regulations. The device has been tested at the manufacturer's site and is delivered in a state safe for operation.

Basic hazards are residual risks that may arise despite proper and safety-conscious use of the device.

DANGER

Risk of death by electrocution!

Body contact with live parts may lead to death or serious injuries.

- Turn off the mains voltage before any works on the device.
- Secure the device against unintentional restart.
- Works on the device may only be executed by service technicians.
- Never operate the device when the mains cable is damaged.

WARNING

Risk of injury due to improper use!

Improper use of the device may lead to injury.

- Always protect the device from extreme heat (excessive sun exposure, close proximity to open fire or heating equipment) during operation and storage.
- Avoid hard impacts that might damage parts of the device.

Staff and qualifications

Permissions Works on/with the device may only be performed by technical staff and/or service technicians (🔑 see "User groups of this document" on page 6).

Note the legally prescribed minimum age for personnel.

Only qualified electricians or technical staff under guidance and supervision of a qualified electrician may perform works on electrical components of the device in accordance with the electrotechnical regulations.

Staff undergoing training or instructions or persons taking part in general vocational training programs may only operate the device under the continuous supervision of technical staff.

Responsibility of the operating company

Condition of the device and/or overall system The operating company shall ensure that all safety and the protective devices are fully functional at all times. This equipment should therefore be checked periodically for functionality and completeness.

This applies to the device as single component as well as the overall system into which the device is integrated as a subcomponent.

Internal measures to avoid hazards The operating company must be aware of the applicable industrial safety regulations and use a risk assessment to determine any additional hazards that may arise from the special working conditions at the device's site of operation. Internal instructions for avoiding the identified hazards should then be drawn up.

Throughout the entire operating period of the device, the operating company must examine and determine whether the supplied operating instructions comply with the current status of applicable regulations.

Staff Regarding the staff authorized or trained by the operating company, the operating company carries the following responsibilities:

- The necessary training and instruction of staff must be guaranteed.
- The powers and responsibilities of staff for installation, operation, maintenance and service must be clearly defined and documented.
- This document must be kept in the immediate surroundings of the device and must be readily accessible to staff.

Working in a safety-conscious manner

Accident prevention and environmental protection In addition to the information in this document, also note the generally applicable statutory and other binding regulations relating to accident prevention and environmental protection.

This may include, for example:

- dealing with hazardous substances
- wearing the necessary and prescribed personal protective clothing
- observance of and compliance with all national and regional industrial safety regulations
- observance of and compliance with all internal working, operating and safety regulations

Content of this document Prior to operation, all staff members assigned to work on the device must have read and understood the relevant parts of this document, especially the chapter “Safety regulations”.

Modifications and alterations

No unauthorized modifications Any unauthorized modifications to the device exclude any liability by the manufacturer for resulting damage and consequences.

Do not perform any modifications, additions and/or alterations to the device without written authorization from point electronic GmbH.

Spare parts and accessories Spare parts and accessories must comply with the technical requirements specified by point electronic GmbH and its suppliers. Whenever original parts are used, compliance is given.

No modifications of the software No modifications to the software must be made or commissioned to third parties. The software may not be deleted, decrypted or decompiled in full or in part.

If changes to the software are necessary, please contact point electronic GmbH.

Maintenance work

Implementation High reliability as well as low maintenance costs of the device may be achieved through careful implementation of the maintenance schedule and regular monitoring over the entire operational lifetime.

Compliance with regulations When performing maintenance work, it is essential to observe the following:

- applicable accident prevention regulations,
- applicable environmental protection regulations and
- safety instructions for maintenance work.

Only perform maintenance work if:

- you are authorized to do so and
- the device is turned off and secured against unintentional restart.

Before getting started Before you begin maintenance work, please make sure that:

- the device is de-energized and disconnected from the power supply and
- the power supply may not be re-established unintentionally as long as the maintenance work lasts.

Works during operation In any case, the operating company or the personnel employed by it must check based on the specific local conditions that the specified work may be performed during operation without endangering anybody.

Continuation next page ...

Electrical equipment Regularly check the electrical equipment of the device. Immediately rectify defects (such as loose cable connections and/or faulty cables).

Damaged or faulty cables may only be replaced by cables that match the requirements stipulated by point electronic GmbH and/or its suppliers!

Check that the surfaces of all electrical equipment parts are dry and free of oil, grease, deposits and corrosion.

Only use voltage-insulated tools!

Do not place any tools or working materials on the conductive surfaces of components.

Device fuses Faulty fuses may only be replaced by fuses that match the requirements (specifications and ratings) stipulated by point electronic GmbH!

Faulty fuses may not be repaired or bypassed, but must be replaced by fuses of the same type.

Cleaning

- Suitable cleaning products** Only use cleaning products approved by point electronic GmbH.
- All cleaning products used by the customer should be checked with point electronic GmbH or the respective supplier to ensure they are compatible with the materials and colors used.
- Unsuitable cleaning agents** Do not clean the device with:
- scratchy, aggressive cleaning agents containing solvents, gasoline, or alcohol,
 - pressured air, high-pressure cleaner or other kinds of cleaning machine.
- Cleaning equipment** Use lint-free cleaning cloths/wipes for cleaning the device.
- To remove stubborn dirt, use lint-free cleaning cloths/wipes moistened with clear water.
- After cleaning** After cleaning, make sure that:
- cables, ports and fittings are free of oil and/or cleaning agents and
 - cables, wires, connectors and electrical components are dry.

Environmental protection

Recyclable materials When performing maintenance work, ensure that reusable materials are recycled.

Disposal Electrical and electronic waste may constitute a hazard to health and environment when disposed off improperly. According to WEEE Directive 2012/19/EU (Waste Electrical and Electronic Equipment Directive), electrical and electronic waste may not be disposed off as general domestic waste but must be handed in at specific collection points or sent back to the manufacturer.

Check with your local environmental protection agency for the prescribed disposal options for commercially used electronic waste.

Only hire waste disposal companies that are approved by the national and regional authorities.



3 System overview

Chapter overview

Purpose This chapter describes the device. It contains information on the use, scope of delivery, equipment, structure and function.

Contents This chapter contains the following information:

› Intended use	23
› Improper use	24
› Scope of delivery.	25
› Optional equipment	26
› System environment requirements	27
› Front panel	28
› Back panel	29
› Touch display	31
› Device labeling	34

Intended use

Purpose The device and software are exclusively intended for the acquisition and management of digital image data.

Operation The device may only be operated in a technically perfect condition and as intended, in compliance with the user manual and under consideration of safety and potential hazards. Malfunctions, in particular those that may impair safety, should be rectified immediately.

If the device shows damages or defects that compromise operational safety, this must immediately be reported to the operations center, and the device must not be put into operation.

Operation and maintenance of the device may only be performed by technical staff and service technicians (🔑 see "User groups of this document" on page 6). The valid safety and accident prevention regulations must be observed.

Compliance with regulations All operating, maintenance, setup and service measures prescribed by the manufacturer must be complied with.

Additional information In addition to this document, the generally applicable, legal and other binding regulations and legislation and environmental protection instructions must be adhered to.

Improper use

Any use not authorized by the manufacturer is not permissible and may lead to injury or property damage.

point electronic GmbH does not accept liability for damage arising from improper use of the device.

In particular, improper use includes:

- operation in explosive environments
- operation in an environment that does not meet the stipulated requirements
 -  see “Operating conditions” on page 60
- operation with opened device housing
- operation with visible damage to the device and/or peripherals (e.g. connected cables, etc.)
- modifications and/or additions that impair the performance
- exchanging components with unauthorized components

Scope of delivery

The following basic equipment is part the scope of delivery:

Component	Item number
DISS 6 Hardware (device)	depends on the type of device delivered
AC power cable	
USB cable (Type-A to Type-B)	
Set of connecting cables for scan and video	
USB stick with: <ul style="list-style-type: none">– Software,– Documentation and– USB device driver	

NOTICE

The complete scope of delivery depends on the respective order.

Optional equipment

The delivered equipment depends on the order details. It includes:

- Multi-Channel Signal Amplifier (MICS)
- Systems for electrical analysis in the SEM

System environment requirements

Minimum requirements The following table contains information on the minimum requirements of the system environment:

Component	Requirement
Computer	IBM-compatible from Core i3 Recommendation: Core i5 <hr/> RAM: 8 GB <hr/> Operating system: from Windows 10 (32/64 Bit) <hr/> Resolution at least 1280×1024 pixel, true color Recommended resolution: 1920×1200 pixel <hr/> At least one free USB 2.0 or USB 3.0 slot Optional: one free LAN slot (RJ-45) for TCP/IP connection
Power connection	At least one socket: – 115/230 V AC – 50/60 Hz – Single phase – Same grounding as the microscope
Periphery	Mouse with scroll wheel

Front panel

Structure The following figure shows the front panel of the device with its components:




Components The following table contains information on the components on the front panel:

No.	Component
1	Touch display <ul style="list-style-type: none"> – displays status and system information for the device – enables various configurations for the device ⓘ see “Touch display” on page 31

Back panel

Structure The following figure shows the back panel of the device with its components:

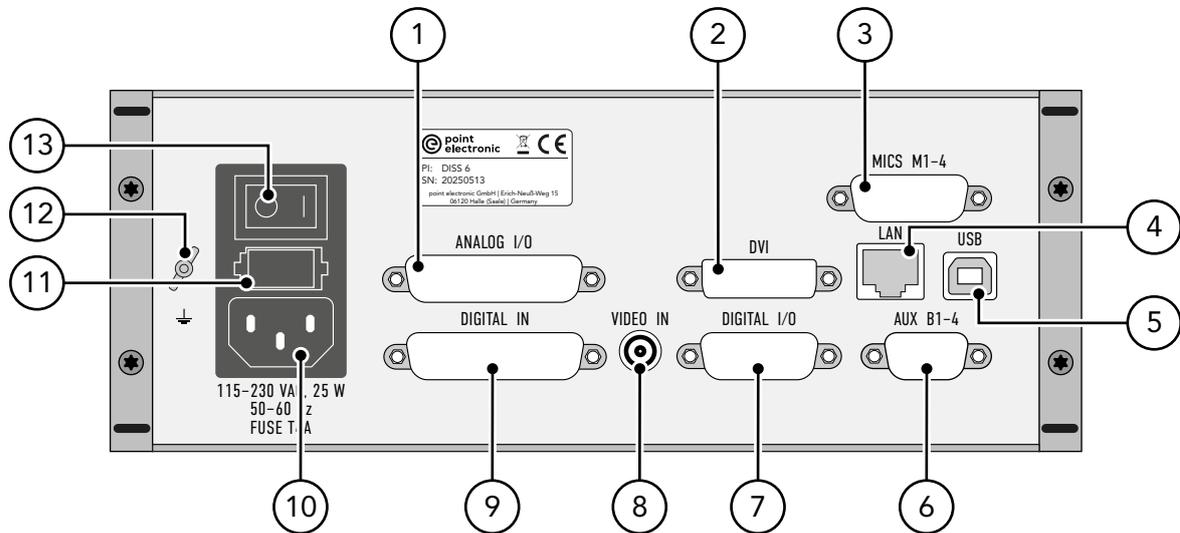

Components The following table contains information on the components on the back panel:

No.	Component
1	<p>ANALOG I/O</p> <ul style="list-style-type: none"> – port for analog inputs and outputs – type: 25 pin, D-Sub, plug (male) <p>Analog outputs</p> <ul style="list-style-type: none"> – amplitude: max. ± 12 V – output offset: ± 2 V <p>Analog inputs</p> <ul style="list-style-type: none"> – max. ± 5 V – ± 0.5 V (with gain = 0 dB) – gain: $-22 \dots +26$ dB – input offset: ± 1.25 V <p> pinout: page 37</p>
2	<p>DVI</p> <ul style="list-style-type: none"> – port – type: DVI-I (dual link), socket (female)
3	<p>MICS M1-4 (optional)</p> <ul style="list-style-type: none"> – input port for Multi-Channel Signal Amplifier (4 channels) – max. ± 5 V – ± 0.5 V (with gain = 1x) – gain: 1 ... 1800x – input offset: ± 1.25 V – type: 15 pin, D-Sub, socket (female) <p> pinout: page 42</p>

Continuation next page ...

No.	Component
4	LAN <ul style="list-style-type: none">– port for TCP/IP control interface– transmission of commands and data– type: RJ-45, socket (female)
5	USB <ul style="list-style-type: none">– port for USB control interface– transmission of commands and data– type: USB type B, socket (female)
6	AUX B1-4 <ul style="list-style-type: none">– port for additional analog inputs (fast)– multiplexed inputs (A or B simultaneous)– max. ± 5 V– ± 0.5 V (with gain = 0 dB)– gain: $-22 \dots +26$ dB– input offset: ± 1.25 V– type: 9 pin, D-Sub, plug (male) <p> pinout: page 41</p>
7	DIGITAL I/O <ul style="list-style-type: none">– port for digital inputs and outputs– input: 3.3/5 V TTL– output: 5 V TTL– type: 15 pin, D-Sub, socket (female) <p> pinout: page 40</p>
8	VIDEO IN <ul style="list-style-type: none">– video input for composite video (PAL)– type: cinch, socket (female)
9	DIGITAL IN <ul style="list-style-type: none">– video counter inputs 1 ... 12– 3.3/5 V TTL– type: 25 pin, D-Sub, socket (female) <p> pinout: page 39</p>
10	Mains voltage connector <ul style="list-style-type: none">– 115 ... 230 V AC, 25 W, 50 ... 60 Hz– type: C14, male
11	Fuses <ul style="list-style-type: none">– 2 pieces– replaceable– type: T4A
12	Grounding connector
13	On/off switch

Touch display

Description The touch display shows status and device information on four switchable pages and allows various configurations for the device.


Status The Status page displays status information about the current scan.

The following figure shows the touch display with the Status page:



Option The Options page shows an overview of the installed hardware options and the size of the pixel map.

The following figure shows the touch display with the Option page:

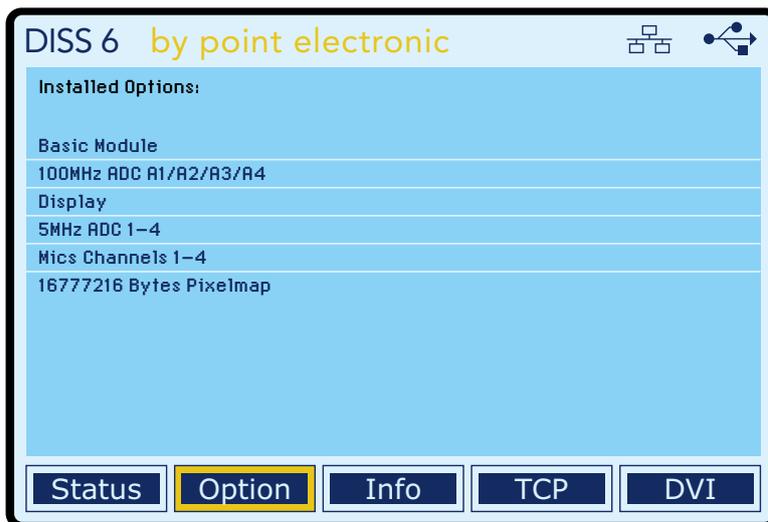

Continuation next page ...


Info The Info page displays information about the installed firmware and current device parameters.

The following figure shows the touch display with the Info page:




TCP The TCP page shows an overview of the current network configuration and allows enabling/disabling of TCP and DHCP.

The following figure shows the touch display with the TCP page:

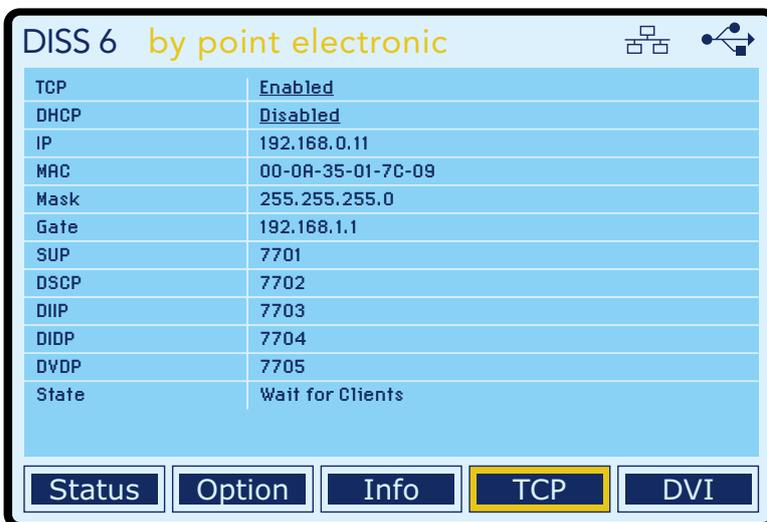

DVI On the page DVI, the DVI output of the device can be configured. This is useful if a monitor is to be connected directly to the device.

The following figure shows the touch display with the DVI page:

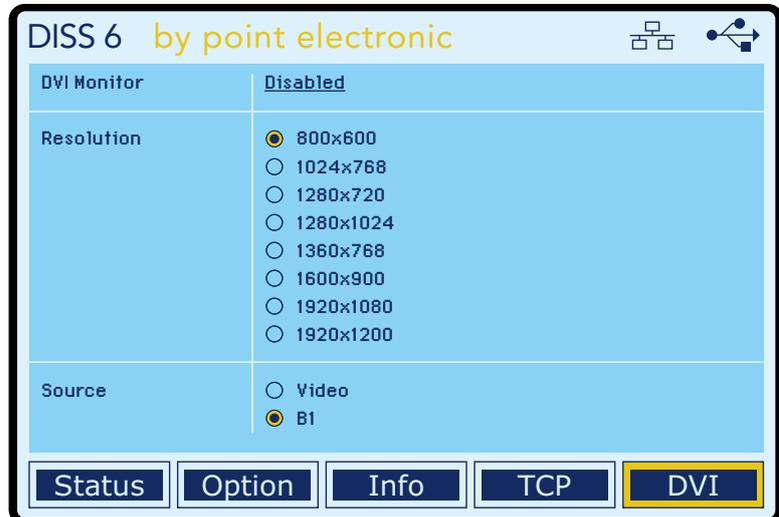


The following options are available for configuring the DVI output:

Option	Description
DVI Monitor	Enabling/disabling the DVI output.
Resolution	Setting the screen resolution.
Source	Setting the input signal source to be displayed on the screen.

Device labeling

Markings on the device The following figure shows the markings on the back of the device:

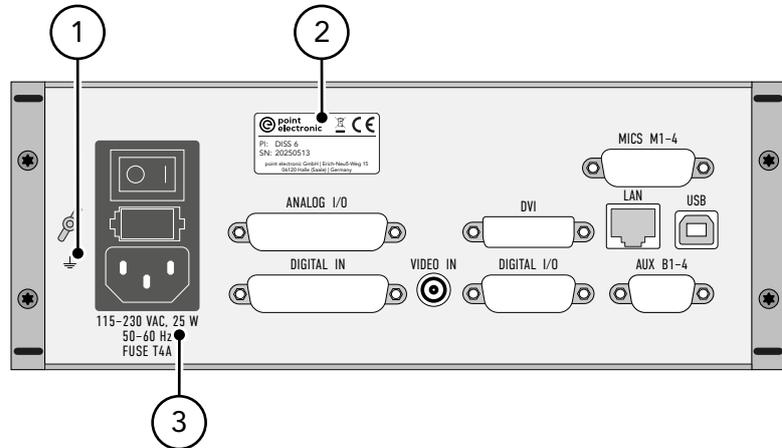


The following table contains information on the markings on the back of the device:

No.	Component
1	Grounding marking
2	Identification plate
3	Information on power supply, power consumption and fuses

Continuation next page ...

Identification plate The following figure shows the identification plate with its components:

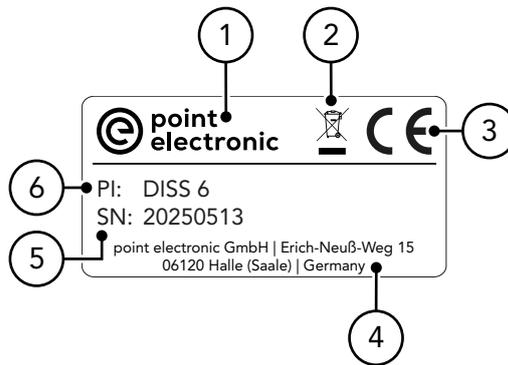

The following table contains information on the components of the identification plate:

No.	Component
1	Manufacturer identification
2	Disposal symbol in accordance with the WEEE directive
3	CE marking
4	Manufacturer address
5	Serial number
6	Device designation



4 Pinouts of the inputs and outputs

Chapter overview

Purpose This chapter describes the pinouts of the particular inputs and outputs.



Contents This chapter contains the following information:

› Analog I/O	37
› Digital in	39
› Digital I/O	40
› AUX B1-4	41
› Mics M1-4	42

Analog I/O

Type 25 pin, D-Sub, plug (male)

Numbering The following table contains information on the numbering of the connections for the respective shape:

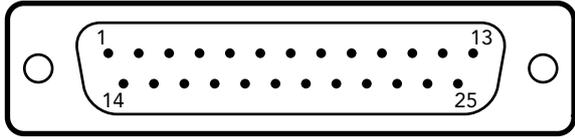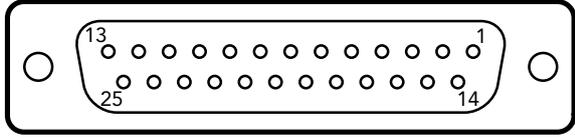Shape	Numbering (view from outside)
Plug	
Socket	

Assignment The following table contains information on the assignment of the connection numbers:

No.	Assignment
1	Scan Y out
2	Scan X out
3	Ext. scan out
4	Ext. blank out
5	Image signal A1+
6	Image signal A1-
7	Image signal A2+
8	Image signal A2-
9	Image signal A3+
10	Image signal A3-
11	Image signal A4+
12	Image signal A4-
13	Not connected
14	Scan Y in
15	Scan X in
16	Ext. scan in
17	Ext. blank in



Continuation next page ...

No.	Assignment
18...24	GND
25	Not connected
Housing	Shield

Digital in

Type 25 pin, D-Sub, socket (female)

Numbering The following table contains information on the numbering of the connections for the respective shape:

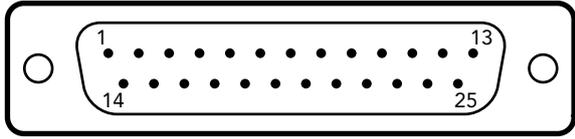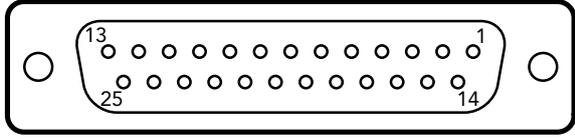Shape	Numbering (view from outside)
Plug	
Socket	



Assignment The following table contains information on the assignment of the connection numbers:

No.	Assignment
1	X 1
2	X 2
3	X 3
4	X 4
5	X 5
6	X 6
7	X 7
8	X 8
9	X 9
10	X 10
11	X 11
12	X 12
13	Not connected
14...25	GND
Housing	Shield

Digital I/O

Type 15 pin, D-Sub, socket (female)

Numbering The following table contains information on the numbering of the connections for the respective shape:

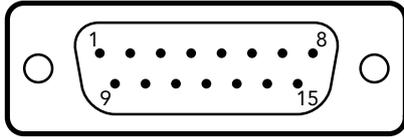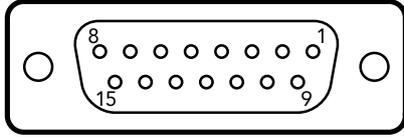Shape	Numbering (view from outside)
Plug	
Socket	



Assignment The following table contains information on the assignment of the connection numbers:

No.	Assignment	Type
1	Pixel clock	output: 5 V TTL
2	Line clock	output: 5 V TTL
3	Frame clock	output: 5 V TTL
4	Hold clock	output: 5 V TTL
5	Ext. scan	output: 5 V TTL
6	Ext. blank	output: 5 V TTL
7	Beam blanker	output: 5 V TTL
8	Lock-In ref out	output: 5 V TTL
9	GND	
10	GND	
11	Pixel sync	input: 3.3/5 V TTL
12	Line sync	input: 3.3/5 V TTL
13	Frame sync	input: 3.3/5 V TTL
14	Hold sync	input: 3.3/5 V TTL
15	Spare I/O	
Housing	Shield	

AUX B1-4

Type 9 pin, D-Sub, plug (male)

Numbering The following table contains information on the numbering of the connections for the respective shape:

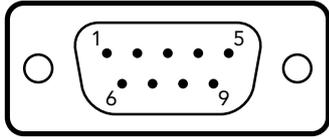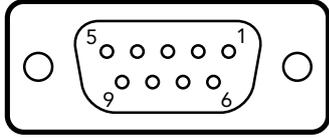Shape	Numbering (view from outside)
Plug	
Socket	



Assignment The following table contains information on the assignment of the connection numbers:

No.	Assignment
1	Image signal B1+
2	Image signal B2+
3	Image signal B3+
4	Image signal B4+
5	GND
6	Image signal B1-
7	Image signal B2-
8	Image signal B3-
9	Image signal B4-
Housing	Shield

Mics M1-4

Type 15 pin, D-Sub, socket (female)

Numbering The following table contains information on the numbering of the connections for the respective shape:

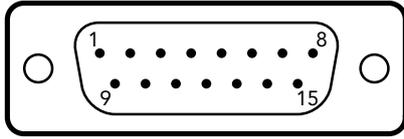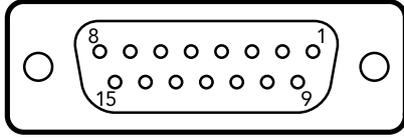Shape	Numbering (view from outside)
Plug	
Shape	

Assignment The following table contains information on the assignment of the connection numbers:

No.	Assignment
1	A+
2	B+
3	C+
4	D+
5	GND
6	+3 V
7	GND
8	+60 V detector BIAS
9	A-
10	B-
11	C-
12	D-
13	GND
14	-3 V
15	GND
Housing	Shield



5 Installation and configuration

Chapter overview

Purpose This chapter contains descriptions and instructions on installation and configuration of the device.

Contents This chapter contains the following information:

- › Configuring the inputs and outputs 44
- › Positioning. 47
- › Connecting to the power supply 48
- › Installing 49
- › Configuring TCP 51

Configuring the inputs and outputs

Description On the carrier board of the device, the inputs and outputs of the device are adjusted to the requirements of the connected SEM.

⚠ DANGER

Risk of death by electrocution!

Body contact with live parts may lead to death or serious injuries as well as the destruction or damaging of components.

- Prior to the configuration of inputs and outputs, please make sure that the device is de-energized and secured against unintentional starting (restart).


Before getting started Complete the following steps before configuring the inputs and outputs of the device:

1. Switch off the device.
2. Remove the mains cable.
3. Remove all connected cables from the device.

Steps Complete the following steps to configure the inputs and outputs of the device:

1. Remove the 4 screws (TORX TX 10) on the back of the device.

 screwdriver TX 10

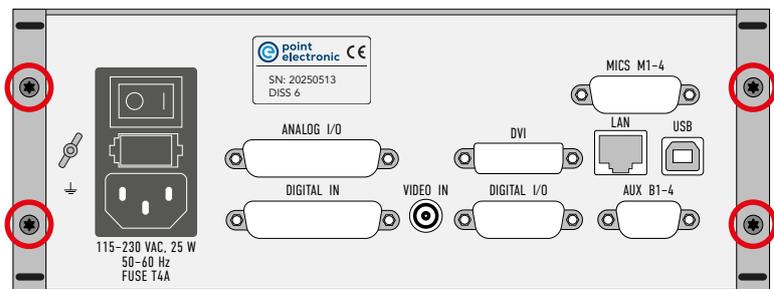


2. Push back the upper casing cover.

Continuation next page ...

... Continuation: Configuring the inputs and outputs

3. Configure the inputs and outputs.
 -  see "Connector strips and assignment" on page 45
 -  see "Configurations for K13 and K14" on page 45
4. Close the casing cover.
5. Attach the 4 screws (TORX TX 10) on the back of the DISS 6 Hardware.
 -  screwdriver TX 10

Connector strips and assignment The following figure shows the distribution of the connector strips and their standard assignment:

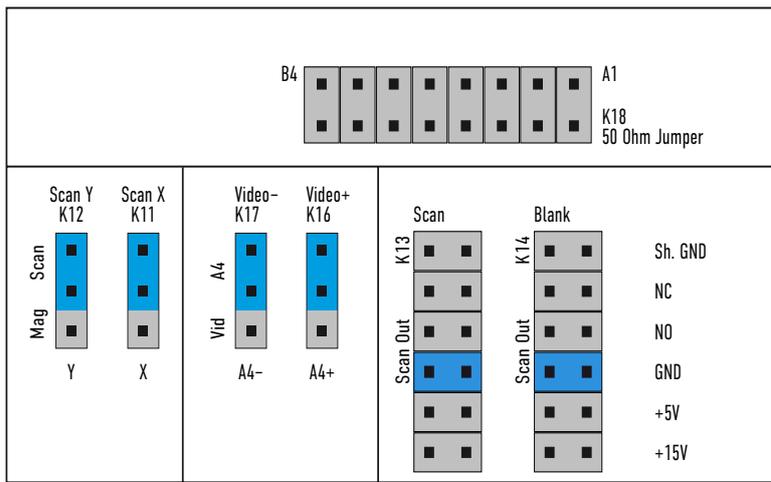

The following table contains information on function and assignment of the connector strips:


Connector strip	Function	Assignment
K18	Terminating resistor for signal inputs A1 ... A4 and B1 ... B4	Jumper closed = terminated with 50 Ω
K11, K12	Type of scan output	<ul style="list-style-type: none"> – Scan (standard) = fixed configured signal level – Mag = signal level adjustable with DAC
K16, K17	Type of signal input A4	<ul style="list-style-type: none"> – A4 (standard) = analog signal – Vid: configuration for composite video (PAL)
K13, K14	Configuration of outputs "ext. Scan" and "ext. Blank"	<ul style="list-style-type: none"> – K13 = signal for ext. Scan – K14 = signal for ext. Blank

Configurations for K13 and K14 The digital outputs "ext. Scan" and "ext. Blank" are potential-free relay outputs destined for switching the SEM to external beam control. They must be adapted to the

Continuation next page ...

requirements of the SEM using the K13 and K14 connector strips.

The following figure shows the possible configurations of the "ext. Scan" and "ext. Blank" outputs on the connector strips K13 and K14:

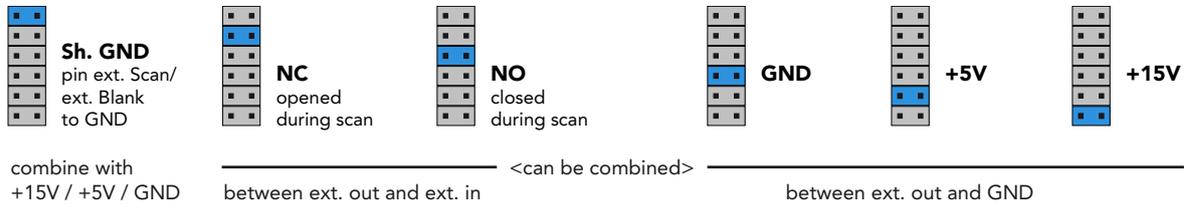

Positioning

Installation site Please make sure that the installation site adheres to the stipulated operating conditions.

 see "Operating conditions" on page 60

Position the device on a level non-slip working surface.

NOTICE

Always ensure accessibility!

When positioning the device, please ensure that on-off switch and all connectors are accessible without obstructions.

Connecting to the power supply

Steps Complete the following steps to connect the device to the power supply:


NOTICE

Mind the device connectors!


When connecting the device, please refer to the description of the unit's backside.

 see "Back panel" on page 29

1. Connect the grounding to the device.

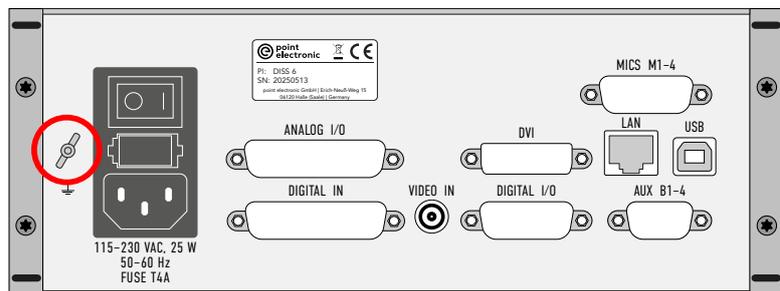

2. Connect the AC power cable to the device.

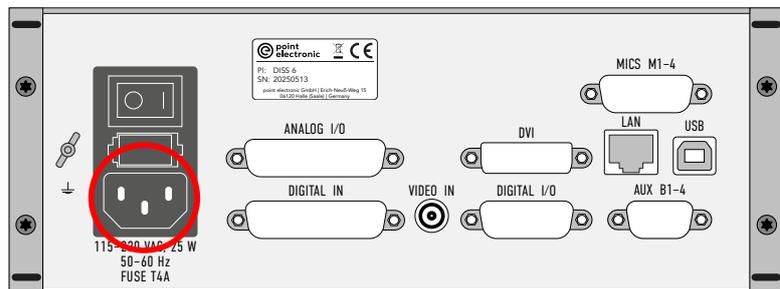

 **Use the AC power cable included in the scope of delivery or an AC power cable that meets the same requirements!**

3. Connect the AC power cable to a power supply.

 **The power supply must comply with the specified requirements!**

 see "Device specifications" on page 61


Installing

Requirements Before installing the device, please make sure that the following requirements are met:

- The inputs and outputs of the device are configured to meet the requirements of the SEM.
 - 📖 see “Configuring the inputs and outputs” on page 44
- The device is connected to a power supply.
 - 📖 see “Connecting to the power supply” on page 48
- The device computer to be connected is equipped with a USB 2.0 or USB 3.0 interface. If this is not the case, a USB plug-in card must be installed. The system requirements must be observed for the driver installation.
 - 📖 see “System environment requirements” on page 27
- The DISS 6 software is installed on the device computer to be connected.

Installing the USB driver Complete the following steps to install the USB driver for the device:

1. Connect the device to a USB port on the device PC using a USB cable.

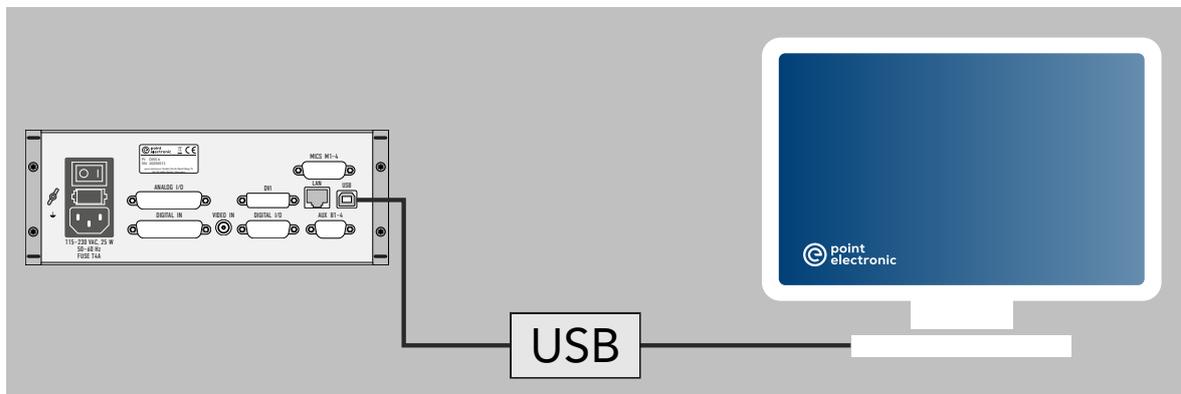

2. Start the device PC.
3. Switch on the device.
 - A dialog window with a prompt to install the driver will appear on the screen.
4. Use the automatic search to find the driver on the device PC.

Continuation next page ...

Integrating into the system Complete the following steps to integrate the device into an existing system:

1. Connect the device to the SEM.

 **Use the connection cables included in the scope of delivery!**

 **Mind the pinouts of the inputs and outputs!**

 see „4 Pinouts of the inputs and outputs“ from page 36

2. Optional: Connect the device to the MICS.
3. Optional: Establish a connection for the camera's synchronization signals.

NOTICE

Perform a functional test!

After complete installation of the DISS 6 Hardware, a functional test may be carried out independent from the SEM.

For this, the X and/or Y deflecting voltage are assigned to the inputs for analog image signals. An image acquisition provides an optical gray wedge for X or Y direction.

Configuring TCP

Description After initial setup, the device may be integrated into the laboratory or microscope network. For the integration, the device's network settings must be configured or adjusted. This is done using the following software:

Setup for ...	Software for setup	Description of the software
EO control	SEM Control Demo	<ul style="list-style-type: none"> – Service Software – Runs on Windows
Imaging	DISS6Test	<ul style="list-style-type: none"> – Can be started directly from a data carrier (e.g., USB stick) without installation

Protocol ports The controllers use different ports for specific network communication protocols.

The default settings for these ports are usually sufficient, unless these ports are already reserved for other communication purposes within the local network.

The following table contains information on the default settings for the particular ports:

Port number	Protocol	Protocol description
7701	SUP	SEM Upgrade Protocol
7702	DSCP	DISS Scan Commissioning Protocol
7703	DIIP	DISS Information Instruction Protocol
7704	DIDP	DISS Image Data Protocol
7705	DVDP	DISS Video Data Protocol

Initial setup via USB For the initial setup, it is recommended to connect the particular device to the device PC via USB. This type of setup is described below.

ATTENTION

Malfunction due to incorrect configuration!

If the TCP configuration is incorrect, the particular device cannot be accessed on the network.

- Before you begin configuration, determine the required settings with a network administrator.

Continuation next page ...

Before getting started Before you begin setting up the TCP connection, ensure that:


- the device is connected to the device PC via USB and switched on and
- the point electronic USB driver is installed on the device PC

Steps Complete the following steps to set up the TCP connection:

1. Start the software "DISS6Test" from the USB stick on the device PC.

i Path to the executable file:
USB Stick > DISS6Test > DISS6Test.exe

2. Open the Misc tab.

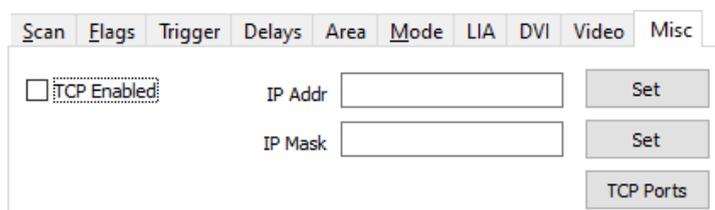

The screenshot shows the 'Misc' tab of the DISS6Test software. The 'TCP Enabled' checkbox is checked. Below it are two input fields: 'IP Addr' and 'IP Mask', each with a 'Set' button to its right. At the bottom right of the tab is a 'TCP Ports' button.

3. Tick the TCP Enabled checkbox to enable TCP.
4. Enter the IP address in the field IP Addr.
⚠ Ensure that the IP address is compatible with the existing settings of the laboratory or microscope network!
5. Click the **Set** button to confirm and save the settings.
6. Enter the IP mask in the field IP Mask.
⚠ Ensure that the IP mask is compatible with the existing settings of the laboratory or microscope network!
7. Click the **Set** button to confirm and save the settings.
8. Optional: Click the **TCP Ports** button to check and/or change the default numbers of the protocol ports.
📖 see "Protocol ports" on page 51
9. Open the Settings menu.
10. Click the menu entry Save Persistent!

Continuation next page ...

11. Connect the device to the laboratory or microscope network.
 -  Only use RJ-45 cable with the specifications CAT.6 and S/FTP (Shielded Foiled Twisted Pair).
12. Restart the device to apply the settings.
13. Open the TCP page on the touch display to check the set IP address.



6 Maintenance

Chapter overview

Purpose This chapter contains instructions on the maintenance of the device.

Contents This chapter contains the following information:

- › Replacing the fuses 55

Replacing the fuses

Position of the fuses The fuses are located at the back of the device.

 TODO: XREF Back panel DISS 6

DANGER

Risk of death by electrocution!

Body contact with live parts may lead to death or serious injuries as well as the destruction or damaging of components.

- Prior to the replacement of the fuses, please make sure that the device is de-energized and secured against unintentional starting (restart).


Before getting started Complete the following steps before replacing the fuses of the device:

1. Switch off the device.
2. Remove the mains cable.
3. Remove all connected cables from the device.

Continuation next page ...

Steps Complete the following steps to replace the fuses of the device:

1. Pinch the locking tongues of the fuse holder and pull the fuse holder from the slot.

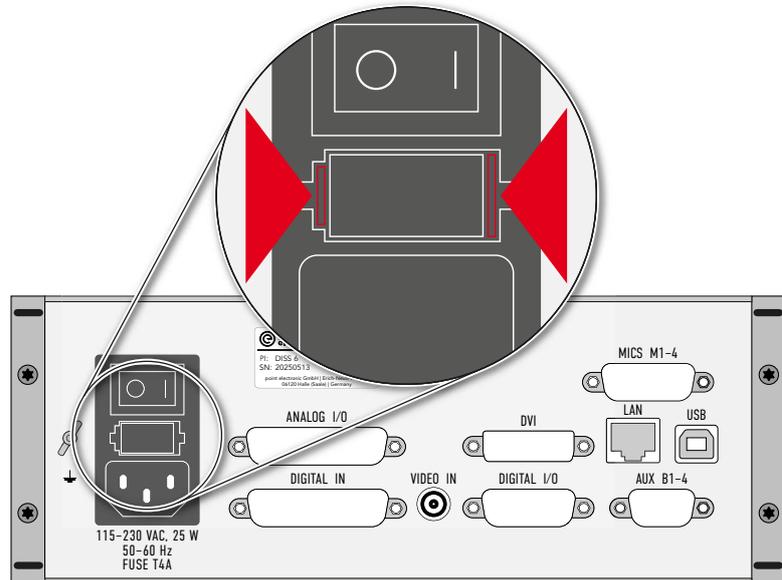

2. Replace the fuse or the fuses.
⚠ Only use type T4A fuses!
3. Push the fuse holder into the slot until the locking tongues snap in.



7 Disposal

Chapter overview

Purpose This chapter contains information on the disposal of the device.

Contents This chapter contains the following information:
› Recycling and taking back of used equipment 58

Recycling and taking back of used equipment

Disposal in accordance with regulations

Electrical and electronic devices may pose a risk to health and the environment if disposed of incorrectly. They cannot therefore be disposed of as domestic waste according to WEEE Directive 2012/19/EU (Waste Electrical and Electronic Equipment Directive). Instead they must be taken to designated collecting points or returned to the manufacturer.

Electrical and electronic devices must undergo specified recycling processes (e.g. with respect to circuit boards) which enable safe, environmentally compatible re-use or separate disposal of different device elements.

The return of used devices is regulated differently in different places. Find out from your local council about the return conditions for commercially used electronic devices.

The device contains no toxic substances requiring separate identification for disposal such as mercury (Hg), cadmium (Cd), lead (Pb) or chrome 6 (e.g. in circuit boards).

Labeling

The following symbol indicates the legal duty to dispose of electronic devices as stipulated:





8 Specifications

Chapter overview

Purpose This chapter lists the specifications of the device.

Contents This chapter contains the following information:

- › Operating conditions 60
- › Device specifications 61

Operating conditions

The following table contains information on the operating conditions of the device:

Site of operation	<ul style="list-style-type: none">– indoor– permanently installed device (mounted in a 19-inch rack)– controlled electromagnetic environment
Ambient air	<ul style="list-style-type: none">– average air pressure (approx. 1 bar)– dry– free of dust
Ambient temperature	5 ... 40°C
Humidity	max. 80% (at 31°C, non-condensing)
Altitude	max. 2,000 m

Device specifications

Electrical safety The following table contains information on electrical safety:

Protection class	I
Protection type	IP20
Overvoltage category	II
Contamination level	1

Power supply The following table contains information on the specifications of the power supply:

Supply voltage	115 ... 230 V AC
Frequency	50 ... 60 Hz
Power consumption	max. 25 W
Fluctuation of the supply voltage	max. $\pm 10\%$

Fuses The following table contains information on the fuses:


Number	2 pc. (replaceable)
Type	T4A
Rating	4 A slow blowing

Dimensions and weight The following table contains information on the dimensions and weights:

Height	90 mm
Width	235 mm
Depth	290 mm
Weight	approx. 3.4 kg

Continuation next page ...

Dimensional drawing The following figure shows the dimensions of the device:

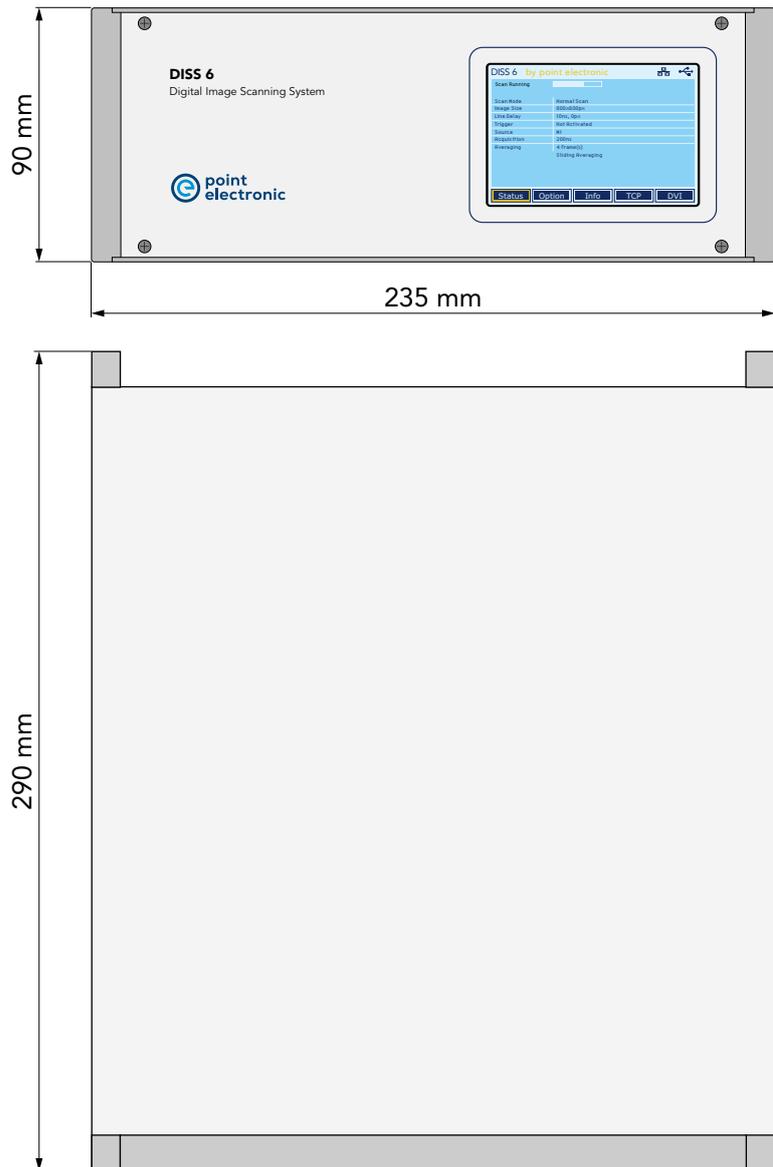

Continuation next page ...

Device housing The following table contains information on the components of the device housing:

Cover plate and base plate	Material: Aluminum Surface: powder-coated RAL 7035 (light gray)
Front panel and rear panel	Material: Aluminum Surface: powder-coated RAL 7035 (light gray)
Frame and profiles	Material: Aluminum Surface: powder-coated RAL 7001 (silver gray)